



Human Elastase ELISA Kit

Vertrieb:

L O X O GmbH Immunbiologie Biochemie, Produkte und Systeme
Postfach 11 30 69215 Dossenheim
Telefon +49 (0) 62 21 - 86 80 23 **FAX** +49 (0) 62 21 - 86 80 255
E-Mail: info@loxo.de **Internet:** www.loxo.de

Assaypro LLC
30 Triad South Drive
St. Charles, MO 63304
T (636) 447-9175
F (636) 447-9475
www.assaypro.com

Hinweis/Note:

Der Packungsbeileger dient nur als erste Information.
Der relevante Packungsbeileger liegt der Ware bei.


The datasheet is only a first information.
The relevant datasheet is included with the product.

For any questions regarding troubleshooting or performing the assay, please contact our support team at support@assaypro.com.


Thank you for choosing Assaypro.

Assay Summary

Add 50 μ l of standard/sample per well.
Incubate 2 hours.

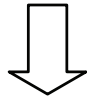

Wash, then add 50 μ l of
biotinylated antibody per well.
Incubate 1 hour.

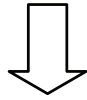

Wash, then add 50 μ l of SP per well.
Incubate 30 minutes.



Wash, then add 50 μ l of
Chromogen Substrate per well.
Incubate 7 minutes.



Add 50 μ l of Stop Solution per well.
Read at 450 nm immediately.

AssayMax Human Neutrophil Elastase ELISA Kit

Catalog No. EE1001-1
Sample Insert/Reference Only

Introduction

Human Neutrophil Elastase (NE), also known as ELA2, is a subfamily of serine proteinase elastases. It consists of 218 amino acid residues with a molecular weight of 23 kDa and two asparagine-linked carbohydrate side chains (1). Neutrophil elastase is stored in neutrophil lysosomes azurophil granules during neutrophil differentiation and is involved in a variety of immune defense reactions and degenerative and inflammatory diseases (2, 3). Upon infection, the activated neutrophils release elastase which then hydrolyzes azurophil granule proteins and extracellular matrix proteins: elastin, collagen, and proteoglycan (4). As a host defense protein, neutrophil elastase degrades bacterial outer membranes and virulence proteins (5, 6). When expressed abnormally, neutrophil elastase causes the development of pulmonary emphysema (7). Neutrophil elastase mutations are associated with cyclic neutropenia and severe congenital neutropenia (8, 9).

Principle of the Assay

The AssayMax Human Neutrophil Elastase ELISA (Enzyme-Linked Immunosorbent Assay) kit is designed for detection of human neutrophil elastase in plasma, serum, urine, saliva, milk, CSF, and cell culture supernatants. This assay employs a quantitative sandwich enzyme immunoassay technique that measures human elastase in less than 4 hours. A polyclonal antibody specific for human elastase has been pre-coated onto a 96-well microplate with removable strips. Neutrophil elastase in standards and samples is sandwiched by the immobilized antibody and the biotinylated polyclonal antibody specific for elastase, which is recognized by a streptavidin-peroxidase conjugate. All unbound material is then washed away and a peroxidase enzyme substrate is added. The color development is stopped and the intensity of the color is measured.

Caution and Warning

- **Prepare all reagents (working diluent buffer, wash buffer, standards, biotinylated antibody, and SP conjugate) as instructed, prior to running the assay.**

- **Prepare all samples prior to running the assay. The dilution factors for the samples are suggested in this protocol. However, the user should determine the optimal dilution factor.**
- **Spin down the SP conjugate vial and the biotinylated antibody vial before opening and using contents.**
- This kit is for research use only.
- The kit should not be used beyond the expiration date.
- The Stop Solution is an acidic solution.

Reagents

- **Human Neutrophil Elastase Microplate:** A 96-well polystyrene microplate (12 strips of 8 wells) coated with a polyclonal antibody against human neutrophil elastase.
- **Sealing Tapes:** Each kit contains 3 precut, pressure sensitive sealing tapes that can be cut to fit the format of the individual assay.
- **Human Neutrophil Elastase Standard:** Human neutrophil elastase in a buffered protein base (20 ng, lyophilized).
- **Biotinylated Neutrophil Elastase Antibody (50x):** A 50-fold concentrated biotinylated polyclonal antibody against neutrophil elastase (140 μ l).
- **EIA Diluent Concentrate (10x):** A 10-fold concentrated buffered protein base (30 ml).
- **Wash Buffer Concentrate (20x):** A 20-fold concentrated buffered surfactant (30 ml, 2 bottles).
- **Streptavidin-Peroxidase Conjugate (SP Conjugate):** A 100-fold concentrate (80 μ l).
- **Chromogen Substrate:** A ready-to-use stabilized peroxidase chromogen substrate tetramethylbenzidine (8 ml).
- **Stop Solution:** A 0.5 N hydrochloric acid to stop the chromogen substrate reaction (12 ml).

Storage Condition

- Store components of the kit at 2-8°C or -20°C upon arrival up to the expiration date.
- Store SP Conjugate and biotinylated antibody at -20°C.
- Store Microplate, Diluent Concentrate (10x), Wash Buffer, Stop Solution, and Chromogen Substrate at 2-8°C.
- Unused microplate wells may be returned to the foil pouch with the desiccant packs and resealed. May be stored for up to 30 days in a vacuum desiccator.
- Diluent (1x) may be stored for up to 30 days at 2-8°C.
- Store standard at 2-8°C before reconstituting with diluent and at -20°C after reconstituting with diluent.

Other Supplies Required

- Microplate reader capable of measuring absorbance at 450 nm.
- Pipettes (1-20 μ l, 20-200 μ l, 200-1000 μ l and multiple channel).
- Deionized or distilled reagent grade water.

Sample Collection, Preparation and Storage

- **Plasma:** Collect plasma using one-tenth volume of 0.1 M sodium citrate as an anticoagulant. Centrifuge samples at 3000 x *g* for 10 minutes. Dilute samples 1:50 into EIA Diluent and assay. The undiluted samples can be stored at -20°C or below for up to 3 months. Avoid repeated freeze-thaw cycles (EDTA or Heparin can also be used as an anticoagulant).
- **Serum:** Samples should be collected into a serum separator tube. After clot formation, centrifuge samples at 3000 x *g* for 10 minutes and remove serum. Dilute samples 1:50 into EIA Diluent and assay. The undiluted samples can be stored at -20°C or below for up to 3 months. Avoid repeated freeze-thaw cycles.
- **Cell Culture Supernatants:** Centrifuge cell culture media at 3000 x *g* for 10 minutes to remove debris. Collect supernatants and assay. Store the remaining samples at -20°C or below. Avoid repeated freeze-thaw cycles.
- **Urine:** Collect urine using sample pot. Centrifuge samples at 800 x *g* for 10 minutes. Dilute urine 1:8 with EIA Diluent and assay. The undiluted samples can be stored at -20°C or below for up to 3 months. Avoid repeated freeze-thaw cycles.
- **Saliva:** Collect saliva using sample tube. Centrifuge samples at 800 x *g* for 10 minutes. Dilute saliva 1:1000 with EIA Diluent and assay. The undiluted samples can be stored at -20°C or below for up to 3 months. Avoid repeated freeze-thaw cycles.
- **Milk:** Collect milk using sample tube. Centrifuge samples at 800 x *g* for 10 minutes. Dilute milk 1:200000 with EIA Diluent or within the range 1:100000 to 1:400000, and assay. The user should determine the optimal dilution factor. The undiluted samples can be stored at -20°C or below for up to 3 months. Avoid repeated freeze-thaw cycles.
- **CSF:** Collect cerebrospinal fluid (CSF) using sample pot. Centrifuge samples at 3000 x *g* for 10 minutes. Dilute samples 1:20 into EIA Diluent and assay. The undiluted samples can be stored at -80°C for up to 3 months. Avoid repeated freeze-thaw cycles.

Reagent Preparation

- Freshly dilute all reagents and bring all reagents to room temperature before use.

- **EIA Diluent Concentrate (10x):** If crystals have formed in the concentrate, mix gently until the crystals have completely dissolved. Dilute the EIA Diluent Concentrate 1:10 with reagent grade water. Store for up to 30 days at 2-8°C.
- **Standard Curve:** Reconstitute the 20 ng of Human Neutrophil Elastase Standard with 2 ml of EIA Diluent to generate a solution of 10 ng/ml. Allow the standard to sit for 10 minutes with gentle agitation prior to making dilutions. Prepare duplicate or triplicate standard points by serially diluting the standard solution (10 ng/ml) 1:2 with EIA Diluent to produce 5, 2.5, 1.25, 0.625, 0.313, and 0.156 ng/ml solutions. EIA Diluent serves as the zero standard (0 ng/ml). Any remaining solution should be frozen at -20°C and used within 30 days.

| Standard Point | Dilution | [Elastase] (ng/ml) |
|----------------|--------------------------------|--------------------|
| P1 | Standard (10 ng/ml) | 10.00 |
| P2 | 1 part P1 + 1 part EIA Diluent | 5.000 |
| P3 | 1 part P2 + 1 part EIA Diluent | 2.500 |
| P4 | 1 part P3 + 1 part EIA Diluent | 1.250 |
| P5 | 1 part P4 + 1 part EIA Diluent | 0.625 |
| P6 | 1 part P5 + 1 part EIA Diluent | 0.313 |
| P7 | 1 part P6 + 1 part EIA Diluent | 0.156 |
| P8 | EIA Diluent | 0.000 |

- **Biotinylated Neutrophil Elastase Antibody (50x):** Spin down the antibody briefly and dilute the desired amount of the antibody 1:50 with EIA Diluent. Any remaining solution should be frozen at -20°C.
- **Wash Buffer Concentrate (20x):** If crystals have formed in the concentrate, mix gently until the crystals have completely dissolved. Dilute the Wash Buffer Concentrate 1:20 with reagent grade water.
- **SP Conjugate (100x):** Spin down the SP Conjugate briefly and dilute the desired amount of the conjugate 1:100 with EIA Diluent. Any remaining solution should be frozen at -20°C.

Assay Procedure

- Prepare all reagents, working standards and samples as instructed. Bring all reagents to room temperature before use. The assay is performed at room temperature (20-25°C).
- Remove excess microplate strips from the plate frame and return them immediately to the foil pouch with desiccants inside. Reseal the pouch securely to minimize exposure to water vapor and store in a vacuum desiccator.

- Add 50 μ l of Human Neutrophil Elastase Standard or sample per well. Cover wells with a sealing tape and incubate for 2 hours. Start the timer after the last sample addition.
- Wash five times with 200 μ l of Wash Buffer manually. Invert the plate each time and decant the contents; hit 4-5 times on absorbent material to completely remove the liquid. If using a machine, wash six times with 300 μ l of Wash Buffer and then invert the plate, decanting the contents; hit 4-5 times on absorbent material to completely remove the liquid.
- Add 50 μ l of Biotinylated Neutrophil Elastase Antibody to each well and incubate for 1 hour.
- Wash the microplate as described above.
- Add 50 μ l of Streptavidin-Peroxidase Conjugate to each well and incubate for 30 minutes. Turn on the microplate reader and set up the program in advance.
- Wash the microplate as described above.
- Add 50 μ l of Chromogen Substrate per well and incubate for about 7 minutes or till the optimal blue color density develops. Gently tap plate to ensure thorough mixing and break the bubbles in the well with pipette tip.
- Add 50 μ l of Stop Solution to each well. The color will change from blue to yellow.
- Read the absorbance on a microplate reader at a wavelength of 450 nm **immediately**. If wavelength correction is available, subtract readings at 570 nm from those at 450 nm to correct optical imperfections. Otherwise, read the plate at 450 nm only. Please note that some unstable black particles may be generated at high concentration points after stopping the reaction for about 10 minutes, which will reduce the readings.


Data Analysis

- Calculate the mean value of the duplicate or triplicate readings for each standard and sample.
- To generate a standard curve, plot the graph using the standard concentrations on the x-axis and the corresponding mean 450 nm absorbance on the y-axis. The best-fit line can be determined by regression analysis using four-parameter or log-log logistic curve-fit.
- Determine the unknown sample concentration from the Standard Curve and multiply the value by the dilution factor.

Standard Curve

- The curve is provided for illustration only. A standard curve should be generated each time the assay is performed.

Human Elastase Standard Curve

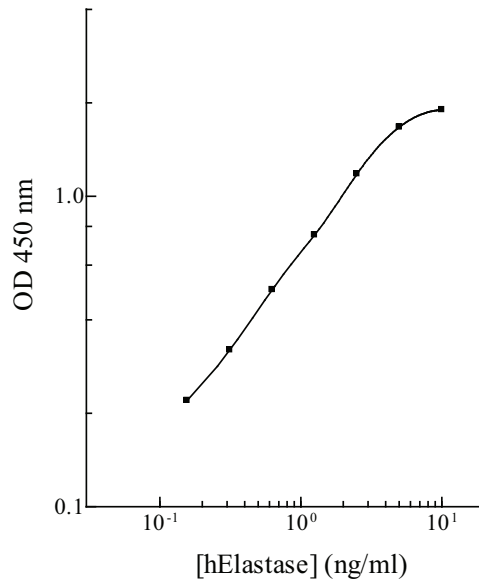

Sensitivity and Specificity

- The minimum detectable dose of elastase is typically ~ 0.15 ng/ml.
- Intra-assay and inter-assay coefficients of variation were 5.0% and 7.1% respectively.

Linearity

| Sample Dilution | Plasma | Serum |
|-----------------|--------|-------|
| 1:25 | 89% | 93% |
| 1:50 | 97% | 100% |
| 1:100 | 101% | 109% |

| Sample Dilution | Saliva |
|-----------------|--------|
| 1:500 | 86% |
| 1:1000 | 97% |
| 1:2000 | 105% |

| Sample Dilution | Urine |
|-----------------|-------|
| 1:4 | 89% |
| 1:8 | 98% |
| 1:16 | 104% |

Recovery

| | |
|-----------------------------|---------------|
| Standard Added Value | 0.5 – 5 ng/ml |
| Recovery % | 84 - 109% |
| Average Recovery % | 96% |

Cross-Reactivity

| Species | % Cross Reactivity |
|----------------|---------------------------|
| Canine | <10% |
| Bovine | None |
| Monkey | <10% |
| Mouse | None |
| Rat | None |
| Swine | None |
| Rabbit | None |
| Human | 100% |

Reference Value

- Normal human elastase plasma levels are <160 ng/ml.

References

- (1) Sinha S *et al.* (1987) *Proc. Natl. Acad. Sci. U.S.A.* 84(8):2228-2232
- (2) Lindmark A *et al.* (1990) *Blood.* 76(11):2374-2380
- (3) Zimmer M *et al.* (1992) *Proc. Natl. Acad. Sci. U.S.A.* 89(17):8215-8219
- (4) Takahashi H *et al.* (1988) *J. Biol. Chem.* 263(29):14739-14747
- (5) Belaaouaj A *et al.* (2000) *Science.* 289(5482):1185-1188
- (6) Weinrauch Y *et al.* (2002) *Nature* 417(6884):91-94
- (7) Guarnieri F *et al.* (2010) *Proc. Natl. Acad. Sci. U.S.A.* 107(23):10661-10666
- (8) Van de Vosse E *et al.* (2011) *Ann Hematol.* 90(2):151-158
- (9) Dale DC *et al.* (2000) *Blood* 96(7):2317-2322

Version 2.4